

FHI 380

FUME CUPBOARD FUNCTION INDICATOR

FHI 380 is an activity indicator to be connected to the controller FHR 320.



Dimensions: 32 x 120 x 18 (6)
(width x height x thickness)

TECHNICAL INFORMATION

Operating voltage	From FHR 320 (6 VDC)
Display	TFT Color Display (135 x 240 pixels)
Switches	Maximum and minimum airflow Alarm acknowledgement Boost and ECO modes (Greenline+ only) Light control (Greenline+ only) Autoclose functions (Greenline+ only)
Cable	2 x 2 x 0,22 mm ² (min)
Enclosure	Polycarbonate IP 44

TOIMINTAKUVAUS

The FHI 350 is a slim operation indicator that can be installed in the fume cupboard glass line or between the fume cupboard and the glass sash.

The control panel is implemented with membrane keys, which allows operation with gloves. The IP44 rated display panel is easy to clean and disinfect.

The switches can be used to forcibly control the control damper to the closed or open position if necessary. When the function is activated, the light on the switch lights up and the display shows the active function.

A air flow alarm occurs if the air flow in the fume cupboard sash opening drops below the set limit. An alarm is indicated by a flashing red light and an alarm sound on the activity indicator (can also be switched off).

CIRCUIT DIAGRAM

Connection to controller FHR 320.

